

STAA SELECTIVE SUPPLEMENT

Vial contents (for 500 ml of medium)

Streptomycin sulphate.....250 mg

Thallium acetate.....25 mg

DIRECTIONS

Aseptically reconstitute the contents of one vial with 5 ml of sterile distilled water. Prepare 500 ml of STAA Agar Base (cat. N° 402079), autoclaved and cooled to 50°C. Add the supplement, mix well and distribute into sterile Petri dishes.

INTENDED USE

For the selective enumeration of *Brochotrix thermosphacta*. For the details of the procedure see the technical sheet of STAA Agar Base (cat. N° 402079)

STORAGE

Store at 2-8° - When stored as directed the supplement remains stable until the expiry date shown on the label. Do not use beyond stated expiry date.

PRECAUTIONS AND WARNING

The supplement should be used only by adequately trained personnel with knowledge of microbiological techniques in the laboratory.

Consult the material safety data sheet before the use.

PACKAGING**4240052 STAA Selective Supplement,****10 vials, each for 500 ml of medium**